

## Iodine-123 Radiochemical Sodium Iodide Solution

### Sodium Iodide (I-123)

**Half-Life: 13.2 hours**

#### IODINE-123 RADIOCHEMICAL SODIUM IODIDE SOLUTION

Iodine-123 is an ideal research tool for Neurology, Oncology and Cardiology.

Researchers prefer using iodine-123 because of the short half-life, clear SPECT images, and because iodine-123 is easy to handle, shield and dispose of.

Nordion has an unparalleled base of technical expertise and production capabilities. State-of-the-art facilities provide a reliable supply of isotopes to customers around the world.


#### PROCESS DATA

##### Nuclear Reaction

Primary



Secondary

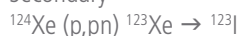

##### Assay

High resolution gamma spectrometry at 159 keV

#### PRODUCT SPECIFICATION

##### Chemical Form

Sodium iodide in 0.1N NaOH

##### Specific Activity

No carrier added

##### Activity Concentration\*

At customer request

##### Radiochemical Purity

≥ 95% iodide

##### Radiopurity\*

≥ 99.8%  $^{123}\text{I}$

≤ 0.2% total other nuclides

\*Activity and radiopurity specifications are at customer's calibration time.

Please note that the radiochemicals listed are not tested for sterility or pyrogenicity. Verification of their suitability for use in humans is the sole responsibility of the purchaser.

#### PACKAGING

- Kontes vial – crimp cap glass V-vial (1 mL, 3 mL, 5 mL)
- Hypovial – crimp cap glass vial with septum (6 mL, 10 mL, 20 mL, 30 mL) (Other vial options are available).

#### AVAILABILITY

Dispensed Monday, Tuesday, Wednesday, Thursday

#### PRODUCT ORIGIN

Vancouver, Canada

#### WORLD CLASS DELIVERY SERVICE

Nordion has a world class distribution system that includes a unique, software-based order management system. Carefully managed transportation logistics ensure products reach their destination on time.



## Iodine-123 Decay Table Half-Life: 13.2 hours

HOURS	0	0.25	0.5	0.75	1	1.25	1.5	1.75	2	2.25	2.5	2.75
0	1.0000	0.9870	0.9741	0.9614	0.9488	0.9365	0.9243	0.9122	0.9003	0.8886	0.8770	0.8655
3	0.8542	0.8431	0.8321	0.8213	0.8105	0.8000	0.7895	0.7792	0.7691	0.7591	0.7492	0.7394
6	0.7297	0.7202	0.7108	0.7016	0.6924	0.6834	0.6745	0.6657	0.6570	0.6484	0.6400	0.6316
9	0.6234	0.6152	0.6072	0.5993	0.5915	0.5838	0.5762	0.5686	0.5612	0.5539	0.5467	0.5396
12	0.5325	0.5256	0.5187	0.5120	0.5053	0.4987	0.4922	0.4858	0.4794	0.4732	0.4670	0.4609
15	0.4549	0.4490	0.4431	0.4373	0.4316	0.4260	0.4204	0.4150	0.4096	0.4042	0.3989	0.3937
18	0.3886	0.3835	0.3785	0.3736	0.3687	0.3639	0.3592	0.3545	0.3499	0.3453	0.3408	0.3363
21	0.3320	0.3276	0.3234	0.3191	0.3150	0.3109	0.3068	0.3028	0.2989	0.2950	0.2911	0.2873
24	0.2836	0.2799	0.2762	0.2726	0.2691	0.2656	0.2621	0.2587	0.2553	0.2520	0.2487	0.2454
27	0.2422	0.2391	0.2360	0.2329	0.2299	0.2269	0.2239	0.2210	0.2181	0.2152	0.2124	0.2097
30	0.2069	0.2042	0.2016	0.1989	0.1964	0.1938	0.1913	0.1888	0.1863	0.1839	0.1815	0.1791
33	0.1768	0.1745	0.1722	0.1699	0.1677	0.1655	0.1634	0.1613	0.1592	0.1571	0.1550	0.1530
36	0.1510	0.1490	0.1471	0.1452	0.1433	0.1414	0.1396	0.1378	0.1360	0.1342	0.1324	0.1307
39	0.1290	0.1273	0.1257	0.1240	0.1224	0.1208	0.1192	0.1177	0.1161	0.1146	0.1131	0.1117
42	0.1102	0.1088	0.1073	0.1059	0.1046	0.1032	0.1019	0.1005	0.0992	0.0979	0.0966	0.0954
45	0.0941	0.0929	0.0917	0.0905	0.0893	0.0882	0.0870	0.0859	0.0848	0.0836	0.0826	0.0815
48	0.0804	0.0794	0.0783	0.0773	0.0736	0.0753	0.0743	0.0734	0.0724	0.0715	0.0705	0.0696
51	0.0687	0.0678	0.0669	0.0660	0.0652	0.0643	0.0635	0.0627	0.0618	0.0610	0.0602	0.0595
54	0.0587	0.0579	0.0572	0.0564	0.0557	0.0550	0.0542	0.0535	0.0528	0.0521	0.0515	0.0508
57	0.0501	0.0495	0.0488	0.0482	0.0479	0.0469	0.0463	0.0457	0.0451	0.0445	0.0440	0.0434

The half-life values stated are as of time of publication. Nordion values may vary.

### CORPORATE HEADQUARTERS

447 March Road  
Ottawa, ON, Canada K2K 1X8  
Tel: +1 613 592 2790  
Fax: +1 613 592 6937

### REGIONAL OFFICE

4004 Wesbrook Mall  
Vancouver, BC, Canada V6T 2A3  
Tel: +1 604 228 8952  
Fax: +1 604 228 5990

### ASIA PACIFIC SALES OFFICES

#### Hong Kong

9/F, 22 Yee Wo Street  
Causeway Bay, Hong Kong  
Tel: +852 2827 8666  
Fax: +852 2827 8302

#### Japan

Room 905, Tokyo Royal Plaza  
1-18-11, Uchi-kanda, Chiyoda-ku  
Tokyo 101-0047, Japan  
Tel: +81 3 5283 6872  
Fax: +81 3 5283 6873

[www.nordion.com](http://www.nordion.com)